

NE-CAT LN₂ FILL STATION

User guide



NECAT LN₂ filling station

Verify all parts of the fill station are connected as shown in the picture. If not, contact your NECAT host.



2. Wear the **PERSONAL PROTECTIVE EQUIPMENT (PPE)**.

PPE (safety glass, face shield, shoe covers, cryogenic gloves and apron) must be worn while operating LN2 filling station.



PURGE

FILL

STOP


“FILL” light should be on. If not, turn the “STOP” button clockwise.

3. NECAT LN₂ control box




4. Fill the Dewar carefully as shown in the picture

Press “FILL” to start filling LN₂. Release the “FILL” button when Dewar is full.

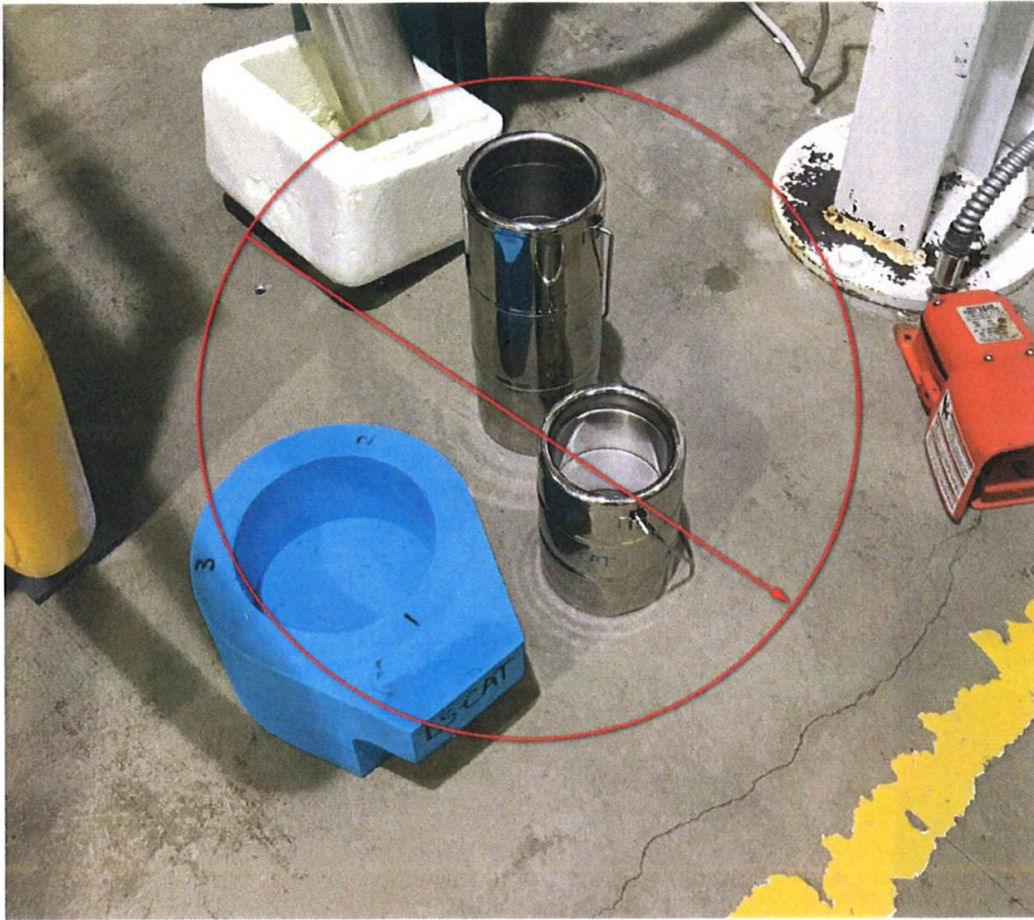

5. STOP. Do not fill small flasks. Use Dewar only

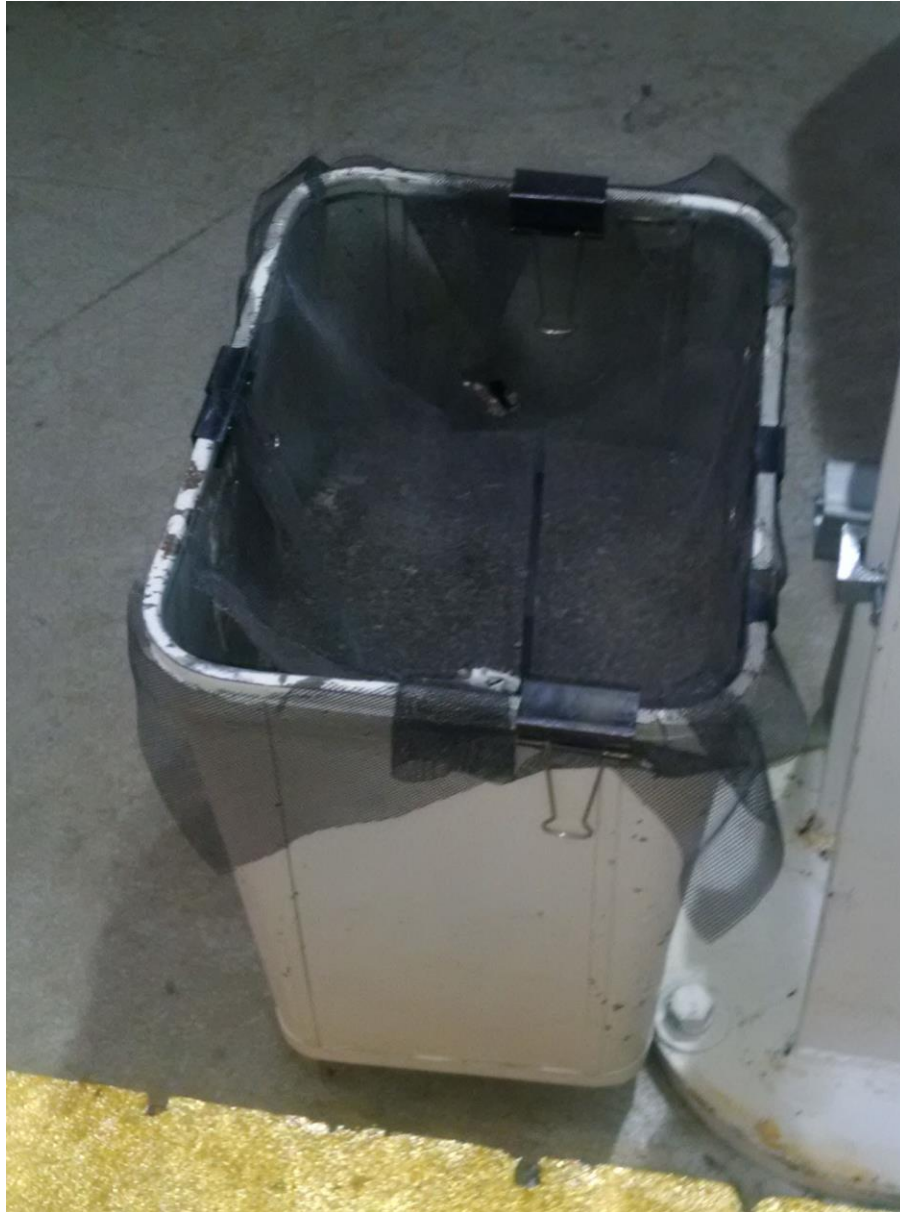

6. Dispose the extra LN₂ in the waste container carefully.

Emergency contacts:

APS Floor coordinator- 2-0101

N. Sukumar 2-0681 (Cell phone 630-328-3532)

Jim Withrow 2-0691 (Cell phone 815-791-4664)

Anthony Lynch 2-0683 (Cell phone 708-620-4190)